

Avant™ buffer pouches

Convenient, pure, fast buffer preparation

Description

Avansta's Avant buffer pouches are perfect for standardizing your electrophoresis and Western blotting applications. Dissolve a pouch of premeasured molecular biology-grade chemicals in 500 ml of deionized water to obtain a ready-to-use 1x buffer without any further preparation. Save precious time in the laboratory and generate reproducible results with Avant buffers.

Ordering Information

R-01038-020	PBS
R-01039-020	TBS
R-01037-020	Tris-Glycine PAGE running buffer
R-01036-020	Tris-Glycine SDS PAGE running buffer



Quick Protocol

1. Open pouch and transfer powder to clean 500 ml container.
2. Add distilled water to a final volume of 500 ml and mix until powder is completely dissolved.
3. Buffer is ready to use.

Tips

1. For best results, use filtered, double-deionized water.
2. There should be no need to adjust pH.
3. For quick identification, each pouch is labeled with a code as follows:

PBS	PBS
TBS	TBS
TG	Tris-Glycine PAGE running buffer
TGS	Tris-Glycine SDS PAGE running buffer

Advansta Corporation

1505 Adams Drive, Suite B1 | Menlo Park, CA 94025

Tel: 650.325.1980 | Fax: 650.325.1904

Email: sales@advansta.com

www.advansta.com

